

Formaldehyde Spill Response Kit Directions

Kit Number: (48625, 48630)

Kit Contents:

- Clear Disposal Bags (Quantity: 4 / Dimensions: 30" x 36")
- FSC 1 Solidifier (11 oz. Bottle)
- Instructions
- Open Back Gown
- Pair of Nitrile Gloves
- Respirator
- Scoop/Scraper (Quantity: 4)
- Twist Ties (Quantity: 4)
- Vented Goggles
- Wiper Pads (Quantity: 3)

Directions:

1. Put on open back gown, respirator, vented goggles and gloves.
*Always wear protective gear and exercise caution during clean up.
2. Sprinkle FSC 1 over spill evenly. The fluid will quickly set; add more FSC-1 as necessary.
3. Remove solidified material with scoop and scraper and place in one of the supplied clear disposal bags.
4. Clean area with a hard surface wipe or disinfectant (For best results use Safetec's[®] SaniZide Plus[®] Germicidal Solution Spray or Germicidal Wipe - sold separately). Wipe up any residue left behind with supplied wiper pads.
3. After procedure is completed remove gloves, vented goggles, respirator and open back gown. Place all contaminated materials in clear disposal bag. Seal bag with supplied twist tie.
4. Dispose of potentially infectious materials in accordance with all local, state and federal regulations.