

Formaldehyde Spill Response Kit Directions

Kit Number: (48625, 48630)

Kit Contents:

- Pair of Nitrile Gloves
- Respirator
- Open Back Gown
- Vented Goggles
- FSC-1 Solidifier (11oz. Bottle)
- Scoop/Scraper (Quantity: 4)
- Wiper Pads (Quantity: 3)
- Clear Disposal Bags (Quantity: 4 / Dimensions: 30" x 36")
- Twist Ties (Quantity: 4)
- Instructions

Directions:

1. Put on open back gown, respirator, vented goggles and gloves.
*Always wear protective gear and exercise caution during clean up.
2. Sprinkle FSC-1 over spill evenly. The fluid will quickly set; add more FSC-1 as necessary.
3. Remove solidified material with scoop and scraper and place in one of the supplied white disposal bags.
4. Clean area with a hard surface wipe or disinfectant (For best results use Safetec's[®] SaniZide Plus[®] surface disinfectant spray or wipe - sold separately). Wipe up any residue left behind with supplied wiper pads.
3. After procedure is completed remove gloves, vented goggles, respirator and open back gown....Place all contaminated materials in white disposal bag. Seal bag with supplied twist tie.
4. Dispose of all waste in accordance with local, state, and federal regulations.