

Instruction Guide

Formaldehyde Spill Response Kit

Reorder # 48625 (Poly Bag) | 48630 (Hard Case)



What's Included:

- 4 Clear Disposal Bags (30" x 36")
- FSC 1 Spill Control (11 oz. Solidifier)
- Instructions
- Pair of Nitrile Gloves
- Open Back Gown
- Respirator
- 4 Scoop/Scrapers
- 4 Twist Ties
- Vented Goggles
- 3 Wiper Pads (11" x 11")



Scan here to view the instruction video!



Step 1: Respirator

Step 1: Put on Formaldehyde Respirator.



Step 2: Gown

Step 2: Put on Open Back Gown.




Step 3: Goggles

Step 3: Put on Vented Goggles.



Step 4: Gloves

Step 4: Put on Nitrile Gloves.

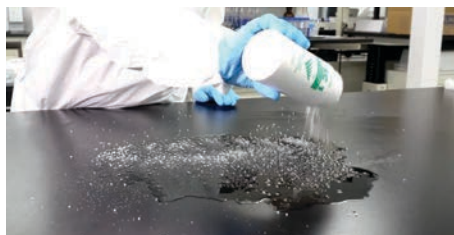

Step 5: FSC 1 Solidifier

Step 5: Sprinkle 11 oz. FSC 1 Solidifier over the spill. Allow liquid to congeal for safer handling and transport.



Step 6: Scoop & Scraper

Step 6: Collect and remove gelled material with Scoop/Scraper.



Step 7: Wiper Pads

Step 7: Wipe up any residue left behind with Wiper Pads.



Step 8: Clear Disposal Bag

Step 8: Place hazardous materials along with Respirator, Protective Gown, Vented Goggles, and Nitrile Gloves into Clear Disposable Bag.



Step 9: Twist Tie

Step 9: Use Twist Tie to securely close Clear Disposable Bag. Dispose according to local, state, and federal regulations.

Rev 25.10