



SPILL CONTROL COLLECTION

FIND YOUR SOLUTION

Meet Safetec 2

Spill Control 3

Red Z® 4-5

Green-Z® 4 & 6

Acid Lock 7

Alky 7

Petro Lock 7

FSC 1 Formalin Spill Control 8

Speedy Cleanz® 8

Zorb Sheets 10

Spill Kits 11

EZ-Cleans Plus® Kit 12

National Standard EZ Cleans® Kit 12

Universal Precaution Compliance Kit 13

EZ-Personal Protection Kit 13

Formaldehyde Spill Kit 14

Hazardous Drugs Spill Kit 14

Multi-Purpose Spill Kit 15

Spill Leader Kit 15

Which Spill Kit is Right for You?..... 16

Spill Kit Extras..... 17

OSHA Compliance 18

5 Steps to Contain & Remove Spills 19

Bloodborne Pathogens Standard 20-21

ANSI Compliance 22

ANSI/ISEA Z308.1-2021 Standard 23-24

ANSI First Aid Kit Checklists 25-26

ANSI First Aid Kit Supplements..... 27-28



Discovering the perfect branding solution has never been easier!

Simplify your search for customization effortlessly with our Private Label Indicator. Spot our symbol in the catalog to identify products eligible for Private Label.

Visit www.privatelabel.safetec.com for more information and to start your journey.




Hello! Welcome to Safetec, where we take pride in offering top-quality infection control, first aid, over-the-counter (OTC), and compliance products. Established in 1992, we have over 30 years of experience in manufacturing. We've made it our mission to uphold the highest quality standards, which is why we strive to be compliant with FDA and EPA regulations and maintain a UL cGMP Certification*. Our diverse product line serves a variety of industries, providing solutions to minimize the risks of cross-contamination and meet safety guidelines, like OSHA and ANSI standards. We are committed to this mission, and we invite you to join us on this journey.

Based in Buffalo, N.Y., we operate three facilities where we oversee every aspect of manufacturing, from concept to product development through distribution. We employ over 120 dedicated professionals in the United States who work tirelessly to manufacture quality products. Whether you're seeking detailed information on our products or need a partner to bring your product to life, we are here to support you every step of the way.

At Safetec, our motto is simple:

One Team, One Goal. Customer Driven.

*Management System Certified by UL with regard to 21 CFR Parts 210/211

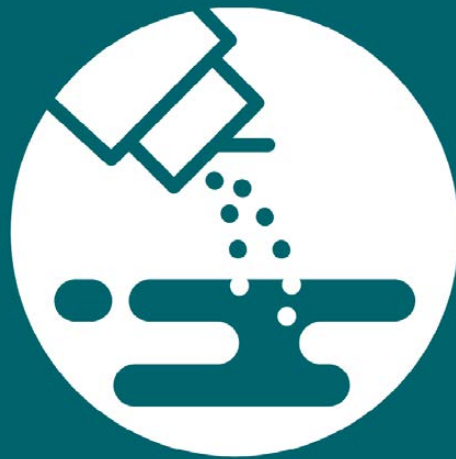

Spill Control

Acid Lock Solidifier	7
Alky Solidifier	7
FSC 1 Formalin Spill Control	8
Green-Z®	4 & 6
Petro Lock Solidifier	7
Red Z®	4-5
Speedy Cleanz®	8
Zorb Sheets	10



Red Z® and Green-Z® Fluid Control Solidifiers



Safetec's Red Z® and Green-Z® solidifiers embody our safety commitment, swiftly containing and solidifying hazardous* spills into a semi-solid (gel) mass. Red Z® handles blood and body fluids, while Green-Z® performs identical functions in addition to chemical spills because there is no chlorine. These solidifiers are user-friendly, eco-conscious, and trusted across industries, and they ensure exceptional performance and regulatory compliance. Available in single & multi-use bottles, single-use pouches, drop-in pacs, and buckets.

- Contains super absorbent polymers (SAPs)
- Eliminates splashing and aerosolization
- Easy-to-use
- Fast-acting
- For potentially infectious blood, body fluid, or chemical spills.***
- Helps comply with DOT regulations†
- Shelf-stable in a cool, dry place**



Deodorizes



Chlorine-Free

**Products are shelf-stable and do not require an expiration date. For the product to remain shelf-stable, it must be kept in a cool, dry place and in its original packaging. If contents are hard/compact, shake the product to loosen before use.

*** Careful consideration should be given to chemical spill cleanup, i.e., room size, temperature, quantity and concentration(s) of the solution(s).

*Dispose of potentially infectious materials in accordance with all local, state and federal regulations.

†Visit our website for a full list of regulation references.



Single & Multi-Use Bottles

Order No.	Size	Absorbency*	Cap	Case Qty.	Case Wt.
41000	1.4 oz.	1,500 cc	Standard	75/case	10 lbs.
41005	2.2 oz.	3,000 cc	Needle Nose	75/case	15 lbs.
41099	7.75 oz.	7,750 cc	Shaker	24/case	14 lbs.
41101	5 oz.	5,000 cc	Shaker	24/case	11 lbs.
41102	11 oz.	11,000 cc	Shaker	12/case	11 lbs.
41103	15 oz.	Up to 18,000 cc	Shaker	12/case	14 lbs.
41105	4.4 oz.	4,400 cc	Needle Nose	24/case	11 lbs.
41107	16 oz.	Up to 22,000 cc	Needle Nose	12/case	14 lbs.

Single-Use & Pour-In Pouches

Order No.	Size	Absorbency*	Bag	Case Qty.	Case Wt.
41109	21 g	787 cc	48 ct.	6/case	15 lbs.
41112	10 g	375 cc	-	500/case	10 lbs.
41117	10 g	375 cc	-	200/case	6 lbs.
41119	21 g	787 cc	-	100/case	6 lbs.
41120	56.7 g	2,126 cc	-	100/case	14 lbs.
41125	21 g	787 cc	-	300/case	16 lbs.
41126	28 g	1,050 cc	-	300/case	21 lbs.
41131	28 g	1,050 cc	-	100/case	8 lbs.

Drop-In Pacs

Order No.	Size	Absorbency*	Bags	Case Qty.	Case Wt.
41108	20 g	750 cc	-	300/case	16 lbs.
41175	40 g	1,500 cc	-	100/case	12 lbs.

Buckets

Order No.	Size	Absorbency*	Product Type	Case Qty.	Case Wt.
41115	3.5 lbs.	56,000 cc	Bucket	1/case	5 lbs.
41116	17.5 lbs.	280,000 cc	Bucket	1/case	22 lbs.
41220	50 lbs.	800,000 cc	Drum	1/case	55 lbs.





Single & Multi-Use Bottles

Order No.	Size	Absorbency*	Cap	Case Qty.	Case Wt.
42003	16 oz.	Up to 22,000 cc	Needle Nose	12/case	15 lbs.
42005	2.2 oz.	3,000 cc	Standard	75/case	16 lbs.
42010	5 oz.	5,000 cc	Shaker	24/case	11 lbs.
42012	15 oz.	Up to 18,000 cc	Shaker	12/case	14 lbs.



Single-Use & Pour-In Pouches

Order No.	Size	Absorbency*	Bags	Case Qty.	Case Wt.
42019	21 g	787 cc	-	100/case	6 lbs.



Drop-In Pacs

Order No.	Size	Absorbency*	Bags	Case Qty.	Case Wt.
42004	10 g	375 cc	100 ct.	10/case	26 lbs.
42008	20 g	750 cc	-	300/case	16 lbs.
42029	4 g	150 cc	500 ct.	10/case	55 lbs.
42030	0.5 g	18.75 cc	500 ct.	3/case	5 lbs.
42031	2 g	75 cc	-	1,000/case	7 lbs.
42175	40 g	1,500 cc	-	100/case	12 lbs.



Buckets

Order No.	Size	Absorbency*	Product Type	Case Qty.	Case Wt.
42220	50 lbs.	800,000 cc	Drum	1/case	55 lbs.

42220 is made to order.



*Data is an approximation of values. Blood and urine solidification are based on the industry standard 0.9% NaCl saline solution.



Acid Lock Solidifier

Rapidly neutralizes and solidifies low pH spills:
Acetic Acid, Hydrochloric Acid, Nitric Acid, and Sulfuric Acid

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
12001	5 oz.	Shaker Top Bottle	-	24/case	11 lbs.
12003	15 oz.	Shaker Top Bottle	-	12/case	14 lbs.



Alky Solidifier

Rapidly neutralizes and solidifies high pH spills:
Ammonium Hydroxide, Sodium Hydroxide, Potassium Hydroxide

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
13001	5 oz.	Shaker Top Bottle	-	24/case	11 lbs.
13003	15 oz.	Shaker Top Bottle	-	12/case	14 lbs.



Petro Lock Solidifier

Rapidly absorbs petroleum-based spills:
Gasoline, Solvents, Crude Oil, and Paint Sludge

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
14101	2 oz.	Shaker Top Bottle	-	24/case	6 lbs.
14103	8 oz.	Shaker Top Bottle	-	12/case	10 lbs.



FSC 1 Formalin Spill Control

Aids in reducing exposure to Formaldehyde Gas by immobilizing formalin and transforming it into an easily removable semi-solid mass. The resulting mass cools rapidly to minimize Formaldehyde off-gassing.

Please Note: Not recommended to be used with 40% formalin or higher.

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
47101	11 oz.	Shaker Top Bottle	-	12/case	11 lbs.

*Looking for a comprehensive solution for Formaldehyde spills?
See page 14 for Safetec's Formaldehyde Spill Response Kit.*



Speedy Cleanz®

Fluid Absorbent

A more economical alternative to our Red Z® Solidifier. Speedy Cleanz® can be applied to vomit, blood, urine as well as other body fluid spills, transforming the liquid into an odorless, semi-solid mass. The non-slippery formula is perfect for use on floors.

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
41100	16 oz.	Shaker Top Bottle	-	12/case	16 lbs.


*Is your facility prepared for a variety of spills?
See Page 15 for the Multi-Purpose Spill Kit and the Spill Leader Kit.*





Zorb Sheets

Blood, Body, & Chemical Fluid Absorbents

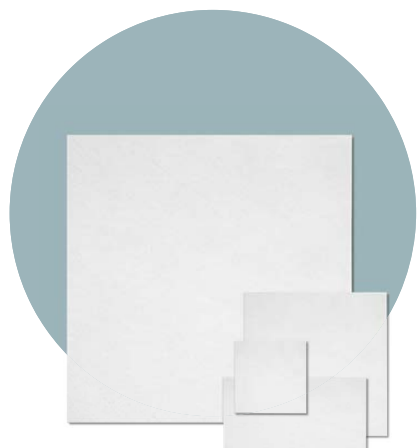

Transfer specimens safely and easily using Safetec's Zorb Sheets! These versatile sheets help reduce compliance risks associated with the handling and shipping of diagnostic specimens, infectious substances, hazardous materials and dangerous goods. The tissue-based material allows fluid to immediately activate the superabsorbent polymers to solidify and transform the contents into a semi-solid (gel) mass. Safetec's Zorb Sheets—your reliable choice for safe transportation that helps comply with DOT, CDC, OSHA, IATA, ICAO, USPS, and UN regulations.†

- Available in multiple sizes
- Fast-acting
- Easy-to-use
- Lightweight
- Space saving
- Prevents leaking

S/S Zorb Sheets

Single-Ply with Standard Absorbency

Order No.	Size	Max. Blood/Urine/Saline Absorbency*	Max. Water Absorbency*	Bags	Case Qty.	Case Wt.
44001	3" x 3"	(Up to 20 cc)	(Up to 131 cc)	1,000 ct.	20/case	26 lbs.
44002	12" x 12"	(Up to 320 cc)	(Up to 2096 cc)	-	500/case	11 lbs.
44003	6" x 6"	(Up to 80 cc)	(Up to 524 cc)	500 ct.	5/case	13 lbs.
44004	3" x 6"	(Up to 40 cc)	(Up to 262 cc)	500 ct.	5/case	7 lbs.

A/S Zorb Sheets

Multi-Ply with Extra Absorbency

Order No.	Size	Max. Blood/Urine/Saline Absorbency*	Max. Water Absorbency*	Bags	Case Qty.	Case Wt.
44016	12" x 12"	(Up to 480 cc)	(Up to 2240 cc)	-	500/case	18 lbs.
44020	2.5" x 4.25"	(Up to 35 cc)	(Up to 165 cc)	100 ct.	50/case	14 lbs.
44035	5" x 6"	(Up to 100 cc)	(Up to 467 cc)	500 ct.	10/case	37 lbs.
44062	4" x 5"	(Up to 67 cc)	(Up to 311 cc)	100 ct.	10/case	5 lbs.
44090	6" x 12"	(Up to 240 cc)	(Up to 1120 cc)	-	1,000/case	18 lbs.
45002	4" x 3"	(Up to 40 cc)	(Up to 187 cc)	1,000 ct.	10/case	31 lbs.
45009	9" x 12"	(Up to 360 cc)	(Up to 1680 cc)	-	500/case	13 lbs.
45010	8" x 9"	(Up to 240 cc)	(Up to 1120 cc)	125 ct.	8/case	17 lbs.

Custom sizes and imprints are available. Call for a quote.

*Data is an approximation of values. Blood and urine solidification are based on the industry standard 0.9% NaCl saline solution.

†Visit our website for a full list of regulation references.



Spill Kits

EZ-Cleans Plus® Kit	12
EZ-Personal Protection Kit	13
Formaldehyde Spill Kit	14
Hazardous Drugs Spill Kit	14
Multi-Purpose Spill Kit	15
National Standard EZ-Cleans® Kit	12
Spill Leader Kit	15
Universal Precaution Compliance Kit	13
Which Spill Kit is Right for You?	16
Spill Kit Extras	17



EZ-Cleans Plus® Kit



The EZ-Cleans Plus® Kit is designed for addressing blood and body fluid spills, and provides essential personal protection items. Its focus on encapsulating and containing potentially infectious spills not only ensures a safer cleanup process but also minimizes the risk of employee exposure to infectious waste and cross-contamination. Each component of the kit is disposable for convenience.

This kit includes Red Z® Solidifier, a quick and efficient solution for solidifying blood and body fluids while eliminating odors.

What's included:

- 10 g Red Z® Solidifier
- Instructions
- Pair of Vinyl Gloves
- Paper Towel
- p.a.w.s.® Antimicrobial Hand Wipe
- Red Biohazard Bag (24" x 24")
- SaniZide Pro 1® Wipe
- Scoop/Scraper
- Twist Tie

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
17120	Plastic Bag	1" x 6.19" x 9.75"	24/case	5 lbs.
17121	Poly Bag	1" x 6.19" x 9.75"	24/case	5 lbs.

17120 does not contain a SaniZide Pro 1® wipe.



Blood and Body Fluid Spill Response



Single-Use

National Standard EZ-Cleans® Kit



Specifically tailored for the cleanup of blood and body fluid spills, the National Standard EZ-Cleans® Kit is a cost-effective alternative for workers who do not require full face and body protection. Compact and easy to use, all kit components are disposable.

This kit includes Red Z® Solidifier, a quick and efficient solution for solidifying blood and body fluids while eliminating odors.

What's included:

- 10 g Red Z® Solidifier
- Instructions
- Odor Mask
- Pair of Vinyl Gloves
- p.a.w.s.® Antimicrobial Hand Wipe
- Red Biohazard Bag (24" x 24")
- SaniZide Pro 1® Wipe
- Scoop/Scraper
- 2 Twist Ties
- White Disposal Bag

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
25000	Refill	1" x 6.19" x 9.75"	24/case	7 lbs.
25005	Hard Case	2.63" x 6.5" x 8.12"	12/case	10 lbs.
25006	Metal Case	2.5" x 5" x 8"	24/case	44 lbs.



Blood and Body Fluid Spill Response



Single-Use



Universal Precaution Compliance Kit



This kit is crucial due to the health risks associated with bloodborne pathogens like HBV, HCV, and HIV. It adheres to OSHA, CDC, and State Health Department standards, providing the necessary personal protection and cleanup tools. Perfect for a multitude of workplaces, including nursing homes, hospitals, schools, correctional facilities, and any business where job-related contact with blood is a potential risk.

This kit includes Red Z[®] Solidifier, a quick and efficient solution for solidifying blood and body fluids while eliminating odors.

What's included:

- 10 g Red Z[®] Solidifier
- Combo Mask/Safety Shield
- Identification Tag
- Instructions
- Pair of Vinyl Gloves
- p.a.w.s.[®] Antimicrobial Hand Wipe
- Protective Apron
- Red Biohazard Bag (24" x 24")
- SaniZide Pro 1[®] Wipe
- Scoop/Scraper
- Twist Tie



Blood and Body Fluid Spill Response



Single-Use

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
17100	Poly Bag	1" x 6.19" x 9.75"	24/case	6 lbs.
17102	Hard Case	2.63" x 6.5" x 8.12"	12/case	10 lbs.

EZ-Personal Protection Kit



This fully disposable kit offers comprehensive exposure protection against potentially infectious blood and body fluids. Designed for pandemic and infection control, this kit comprises a well-organized selection of personal protective equipment (PPE) to address potential exposure to infectious materials. Compact and easily storable, this kit can be safely disposed of after use, ensuring the highest safety and hygiene standards.

This kit includes Red Z[®] Solidifier, a quick and efficient solution for solidifying blood and body fluids while eliminating odors.

What's included:

- Combo Mask/Safety Shield
- Instructions
- Pair of Vinyl Gloves
- p.a.w.s.[®] Antimicrobial Hand Wipe
- Protective Gown
- Red Biohazard Bag (24" x 24")
- Twist Tie



Blood and Body Fluid Spill Response



Single-Use

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
17606	Plastic Bag	1" x 6.19" x 9.75"	24/case	10 lbs.



Formaldehyde Spill Response Kit



A convenient emergency response kit for medical personnel using formalin or other formaldehyde-containing substances. It is ideal for containing small-volume spills.

This kit includes FSC 1, a formalin spill control solidifier that aids in reducing the exposure to formaldehyde gas by immobilizing formalin and transforming it to an easily removable semi-solid mass.

Please Note: Not recommended to be used with 40% formalin or higher.

What's included:

- 4 Clear Disposal Bags (30" x 36")
- FSC 1 Spill Control (11 oz. Solidifier)
- Instructions
- Open Back Gown
- Pair of Nitrile Gloves
- Respirator
- 4 Scoop/Scrapers
- 4 Twist Ties
- Vented Goggles
- 3 Wiper Pads (11" x 11")



Chemical Spill Response



Single-Use

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
48625	Poly Bag	11.75" x 11.75" x 4.88"	1/case	4 lbs.
48630	Hard Case	14" x 12.625" x 5.375"	1/case	6 lbs.

Need to stock up on formaldehyde solidifier? See page 8 for FSC 1 Spill Control.

Hazardous Drugs Spill Kit



In environments where meticulous precautions are paramount like compounding pharmacies and healthcare facilities handling Hazardous Drugs (HDs), such as cytotoxic drugs, Safetec's Hazardous Drugs Spill Kit serves as the primary defense against unexpected spills and exposures.

Any facility that handles HDs must have USP <800> compliant spill kits on-site. This requirement is part of USP General Chapter 800 and is an effort to protect workers and patients in the event of accidental spills of HDs.

What's included:

- 21 g Green-Z® Fluid Solidifier
- Caution Sign
- Chemotherapy Gown
- Full-Face Safety Shield
- Identification Tag
- Instruction Card
- Nitrile Chemotherapy Gloves
- Scoop/Scraper
- Shoe Covers, Non-Slip
- Twist Tie
- 2 Wiper Pads (11" x 11")
- Yellow Chemotherapy Waste Bag (20" x 13" x 39")



Chemical Spill Response



Single-Use

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
49725	Plastic Bag	2" x 8" x 14"	12/case	11 lbs.



Multi-Purpose Spill Kit



This comprehensive kit is your go-to solution for quick and efficient response to a variety of spills. Whether it's chemicals, body fluids, or other hazardous substances, this kit equips you with the essential tools for safe spill cleanup. Compact and easily accessible, it's ideal for a range of environments, from laboratories to workplaces.

Meets requirements for compliance with DOT 1713 hazardous discharge; EPA 623.30 immediate action and 263.31 discharge cleanup.

What's included:

- Acid Lock Solidifier (15 oz. Bottle)
- Alky Solidifier (15 oz. Bottle)
- Green-Z® Solidifier (15 oz. Bottle)
- Petro Lock Absorbent (8 oz. Solidifier)
- Caution Tape Roll
- 3 Clear Disposal Bags (30" x 37")
- D.O.T. Incident Report Tag
- Drop Cloth
- Instructions
- Pair of Nitrile Gloves
- Respirator
- 3 Scoop/Scrapers
- 3 Twist Ties
- Vented Goggles
- Zammie® Sewer Stopper (15" x 21")



Universal Spill Response



Multi-Use

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
15201	Poly Bag	12" x 17.75" x 4.25"	1/case	11 lbs.

Need to refill your solidifiers?
See Page 3 for all Spill Control products.



Spill Leader Kit



This wall-mounted spill kit is designed for emergency response and hazardous lab spills. The solid solutions neutralize, stabilize, and solidify a variety of aqueous spills wherever they occur.

Meets requirements for compliance with DOT 1713 hazardous discharge; EPA 623.30 immediate action and 263.31 discharge cleanup.

What's included:

- Adhesive Wall Mount
- Acid Lock Solidifier (5 oz. Bottle)
- Alky Solidifier (5 oz. Bottle)
- Green-Z® Solidifier (5 oz. Bottle)
- Petro Lock Absorbent (2 oz. Solidifier)
- Red Z® Solidifier (5 oz. Bottle)
- D.O.T. Incident Report Tag
- 2 Hazardous Waste D.O.T Labels
- 3 pairs of Nitrile Gloves
- 2 Non-Hazardous Waste D.O.T. Labels
- 6 Scoop/Scrapers
- 6 Zip Lock Bags (9" x 12")



Universal Spill Response



Multi-Use

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
15205	Poly Bag	14.125" x 3.125" x 20.5"	1/case	5 lbs.

Need to refill your solidifiers?
See Page 3 for all Spill Control products.



Which Spill Kit is Right for You?



Did You Know:

If you have a first aid kit, you need to have a spill kit in order to be OSHA compliant.

Our Universal Precaution Compliance Kit meets OSHA Compliance 1910.1030 Bloodborne Pathogens Standard.

	EZ-Cleans Plus® Kit	National Standard EZ-Cleans® Kit	Universal Precaution Compliance Kit*	EZ-Personal Protection Kit	Formaldehyde Spill Response Kit	Hazardous Drugs Spill Kit	Multi-Purpose Spill Kit	Spill Leader Kit
Cleaning Blood & Body Fluid Spills	✓	✓	✓	✓	-	-	✓	✓
Cleaning Chemical Spills	-	-	-	-	✓	✓	✓	✓
Single-Use	✓	✓	✓	✓	✓	✓	-	-
Multi-Use	-	-	-	-	-	-	✓	✓
Solidifier Included	✓	✓	✓	-	✓	✓	✓	✓
Surface Disinfectant Included	✓	✓	✓	-	-	-	-	-
PPE Included	✓	✓	✓	✓	✓	✓	✓	✓
Portable	✓	✓	✓	✓	✓	✓	✓	-
Stationary	✓	✓	✓	✓	✓	✓	✓	✓

Spill Kit Extras

Elevate your spill control preparedness by having extra items on hand.



Clean Up Essentials

Ensure you are ready for unexpected accidents by maintaining a surplus of essential cleanup supplies like these.


Order No.	Description	Dimensions	Case Qty.	Case Wt.
17170	Scoop/Scraper	-	300/case	8 lbs.
2110122	Red Biohazard Bag	24" x 24"	500/case	16 lbs.
2110105	Red Biohazard Bag	33" x 39"	150/case	12 lbs.
2110014	Poly-Lined Paper Bag (Vomit)	3.25" x 4.75" x 9"	1,000/case	36.5 lbs.



Clean and Disinfect

After taking care of spills, cleaning and disinfecting contaminated areas and maintaining personal hygiene are important.

Order No.	Description	Size	Case Qty.	Case Wt.
13377	Instant Hand Sanitizer 0.9 g Pouch	144 ct.	12/case	7 lbs.
34400B	p.a.w.s.® Antimicrobial Hand Wipe	Bulk	5,000/case	42 lbs.
34810	SaniZide Plus® Germicidal Solution Spray Bottle	32 oz.	6/case	15 lbs.
35910	SaniZide Pro 1® Surface Disinfectant Spray Bottle	32 oz.	6/case	14 lbs.
34452	SaniWash® Antimicrobial Hand Wash	16 oz.	12/case	16 lbs.

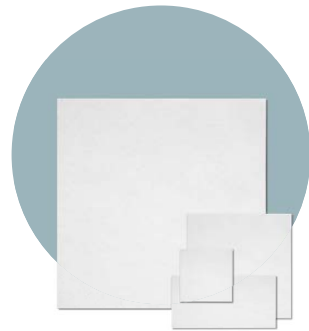

Zorb Sheets

You never know when you may need additional absorption to control a spill. Stock up on Zorb Sheets to be prepared for the unexpected.

Order No.	Description	Bags	Case Qty.	Case Wt.
44001	S/S Zorb Sheets, Standard Absorbency, 3" x 3"	1,000 ct.	20/case	26 lbs.
44002	S/S Zorb Sheets, Standard Absorbency, 12" x 12"	-	500/case	11 lbs.
44003	S/S Zorb Sheets, Standard Absorbency, 6" x 6"	500 ct.	5/case	13 lbs.
44004	S/S Zorb Sheets, Standard Absorbency, 3" x 6"	500 ct.	5/case	7 lbs.

For other available sizes, see Page 10.



Fluid Solidifiers

You can't always anticipate the size of a spill; make sure to have extra solidifier in case your spill is larger than your kit's solidifying abilities.

Order No.	Description	Size	Case Qty.	Case Wt.
41103	Red Z® Shaker Top Bottle	15 oz.	12/case	15 lbs.
41116	Red Z® Bucket	17.5 lbs.	1/case	21 lbs.
41220	Red Z® Drum	50 lbs.	1/case	55 lbs.
42012	Green-Z® Shaker Top Bottle	15 oz.	12/case	15 lbs.
42200	Green-Z® Drum	50 lbs.	1/case	55 lbs.
41100	Speedy Cleanz® Shaker Top Bottle	16 oz.	12/case	16 lbs.
47101	FSC I Formalin Spill Control Shaker Top Bottle	11 oz.	12/case	11 lbs.
12003	Acid Lock Shaker Top Bottle	15 oz.	12/case	14 lbs.
13003	Alky Shaker Top Bottle	15 oz.	12/case	14 lbs.
14103	Petro Lock Shaker Top Bottle	8 oz.	12/case	10 lbs.

42200 is Made to Order.

For other available sizes, see Pages 5-8.



The Occupational Safety and Health Administration's (OSHA) role is to assure America's workers' safety and health by setting and enforcing standards, providing training, outreach, and education, establishing partnerships, and encouraging continual process improvement in workplace safety and health. OSHA sets standards, or rules, that describe the methods that employers must use to protect their employees from hazards.

At Safetec of America, we can offer you the resources and products needed to keep your workplace safe and OSHA-compliant.

5 Steps to Contain & Remove Spills	19
Bloodborne Pathogens Standard	20-21
EZ-Personal Protection Kit	13 & 21
EZ-Cleans® Plus Kit	12 & 21
National Standard EZ-Cleans® Kit	12 & 21
Universal Precaution Compliance Kit ...	13 & 20

5 Steps to Efficiently Contain and Remove Bloodborne Pathogen Spills



29 CFR 1910.1030 – Bloodborne Pathogens

Universal Precautions is an approach to infection control. According to the concept of Universal Precautions, all human blood and certain human body fluids are treated as if known to be infectious for HIV, HBV, and other bloodborne pathogens.

Step 1: Prevent



To prevent direct contact with the spill during the cleanup process and minimize the risk of infection, disposable protective clothing must be worn. This includes gloves, combo mask/safety shield, and a protective apron.

Step 4: Dispose



Discard the germicidal wipe, gloves, safety shield and apron into the **red biohazard bag** and dispose of it according to local, state, and federal regulations.

Step 2: Contain & Remove



To efficiently contain and remove the spill, a **solidifying agent** (Safetec Red Z[®]) and a **scoop/scraper** are to be used. The spill contents are to be disposed of into a **red biohazard bag**.

Step 5: Sanitize



Use an **antimicrobial hand wipe** (p.a.w.s.[®]) to sanitize hands, then use **soap** (SaniWash[®]) and water once the spill is contained and cleaned.

Step 3: Disinfect



To properly disinfect the contaminated surface area, use a **surface disinfectant wipe** (SaniZide Pro 1[®]). This step is critical for preventing further contamination.

Helpful Tips to Remain OSHA Compliant

- Think safety culture – not just compliance
- Compliance is the key to prevention
- Do not guess information or make assumptions
- Plan and prepare for an inspection ahead of time

For more information on OSHA 1910.1030 compliance, visit our website.

Protect against exposure to bloodborne pathogens.



29 CFR 1910.1030 – Bloodborne Pathogens

Universal Precautions is an approach to infection control. According to the concept of Universal Precautions, all human blood and certain human body fluids are treated as if known to be infectious for HIV, HBV, and other bloodborne pathogens.

Following OSHA 29 CFR 1910.1030 is vital for protecting employees from serious health risks, ensuring legal compliance, fostering a safe work environment, and minimizing financial and legal liabilities for employers.

Under this compliance standard, when there is occupational exposure to blood and body fluids, employers are required to provide appropriate, readily accessible personal protective equipment (PPE) at no cost to employees. This PPE must prevent blood or other infectious materials from reaching employees' skin, clothes, eyes, mouth, or mucous membranes during normal use. Contaminated work surfaces must be decontaminated with an appropriate disinfectant after a spill cleanup. Additionally, a means to sanitize hands and other skin must be available, such as antiseptic wipes and soap.

Adhering to these standards not only promotes workplace safety but also helps prevent the transmission of infectious diseases.

With just one kit, the Universal Precaution Kit can help you remain compliant with the OSHA 29 CFR 1910.1030 Bloodborne Pathogens Standard.

Safetec can help!

Universal Precaution Compliance Kit

This kit is crucial due to the health risks associated with bloodborne pathogens like HBV, HCV, and HIV. It adheres to OSHA, CDC, and State Health Department standards, providing the necessary personal protection and cleanup tools.

Suitable for various workplaces, such as nursing homes, hospitals, schools, correctional facilities, and any business where job-related contact with blood is a potential risk.

This kit includes Red Z[®] Solidifier, a quick and efficient solution for solidifying blood and body fluids while eliminating odors.

What's included:

- 10 g Red Z[®] Solidifier
- Combo Mask/Safety Shield
- Identification Tag
- Instructions
- Pair of Vinyl Gloves
- p.a.w.s.[®] Antimicrobial Hand Wipe
- Protective Apron
- Red Biohazard Bag (24" x 24")
- SaniZide Pro 1[®] Wipe
- Scoop/Scraper
- Twist Tie



Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
17100	Poly Bag	1" x 6.19" x 9.75"	24/case	6 lbs.
17102	Hard Case	2.63" x 6.5" x 8.12"	12/case	10 lbs.

Make Any Safetec Kit OSHA Compliant with

EZ-Personal Protection Kit

Combine this PPE bundle with any regular blood and body fluid kit to elevate it to OSHA Bloodborne Pathogen Compliance Standards. This portable, compact kit is perfect for having multiple on hand to handle the unexpected.



What's included:

- Combo Mask/Safety Shield
- Instructions
- Pair of Vinyl Gloves
- p.a.w.s.® Antimicrobial Hand Wipe
- Protective Gown
- Red Biohazard Bag (24" x 24")
- Twist Tie

Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
17606	Plastic Bag	1" x 6.19" x 9.75"	24/case	10 lbs.

Additional Blood and Body Fluid Kits

EZ-Cleans Plus® Kit

What's included:

- 10 g Red Z® Solidifier
- Instructions
- Pair of Vinyl Gloves
- Paper Towel
- p.a.w.s.® Antimicrobial Hand Wipe
- Red Biohazard Bag (24" x 24")
- SaniZide Pro 1® Wipe
- Scoop/Scraper
- 2 Twist Ties



Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
17120	Plastic Bag	1" x 6.19" x 9.75"	24/case	5 lbs.
17121	Poly Bag	1" x 6.19" x 9.75"	24/case	5 lbs.

17120 does not contain a SaniZide Pro 1® wipe.

National Standard EZ-Cleans® Kit

What's included:

- 10 g Red Z® Solidifier
- Instructions
- Odor Mask
- Pair of Vinyl Gloves
- p.a.w.s.® Antimicrobial Hand Wipe
- Red Biohazard Bag (24" x 24")
- SaniZide Pro 1® Wipe
- Scoop/Scraper
- Twist Tie
- White Disposal Bag



Order No.	Container	Container Dimensions	Case Qty.	Case Wt.
25000	Refill	1" x 6.19" x 9.75"	24/case	7 lbs.
25005	Hard Case	2.63" x 6.5" x 8.12"	12/case	10 lbs.
25006	Metal Case	2.5" x 5" x 8"	24/case	44 lbs.



ANSI is the American National Standards Institute and sets guidelines to ensure products and services in the U.S. are safe, reliable, and of high quality. ANSI doesn't make the standards themselves but coordinates the efforts of different groups and companies to develop them. The standards help with consistency and ensure everything works well together, promoting safety and efficiency in various fields.

At Safetec, we manufacture a variety of products that comply with ANSI standards to ensure workplace safety. Our antibiotic ointments, antiseptic, burn treatment, and hand sanitizer lines meet the requirements set by ANSI.

ANSI/ISEA Z308.1-2021 Standard	23-24
ANSI First Aid Kit Checklists	25-26
ANSI First Aid Kit Supplements	27-28

ANSI / ISEA Z308.1-2021 Standard

First Aid Kit Classifications

A

CLASS A KITS

Intended to provide basic range of products to aid most common workplace injuries such as minor cuts, abrasions, and sprains.

B

CLASS B KITS

Intended to provide broader range and quantity of supplies to aid injuries in more complex or high-risk workplace environments.

OSHA 29 CFR 1910.151(b)

OSHA's 29 Code of Federal Regulations (CFR) 1910.151(b) requires that "Adequate first aid supplies shall be readily available." This rule applies to treatment of minor injuries that occur in the workplace. OSHA refers employers to ANSI Z308.1-2021 for guidance on minimum requirements for first aid kits.

ANSI / ISEA Z308.1-2021

American National Standard Minimum Requirements for Workplace First Aid Kits and Supplies Z308.1-2021 establishes minimum performance requirements for first aid kits and their supplies that are intended for use in various work environments. Classification of first aid kits, designating the assortment of items and quantity of each item, is based on the anticipated number of users intended to be served by each first aid kit, as well as the complexity of the work environment and level of hazards.

Safetec® First Aid Supplies

(To help with ANSI / ISEA Compliance Z308.1-2021)

Section 6.4 | Antibiotic Application

Each antibiotic application shall meet the applicable requirements as defined in the current edition of USP/NF. Each application shall be packaged in individual-use, non-reusable applications containing at least 0.14 fl. oz. (0.5 g).

Order No.	Size	Description	Carton	Case Qty.	Case Wt.
53210	0.9 g	Triple Antibiotic Ointment	144 ct.	12/case	7 lbs.
53204	0.9 g	Triple Antibiotic Ointment	Bulk	2,000/case	7 lbs.
53610	0.9 g	Single Antibiotic with Neomycin	144 ct.	12/case	7 lbs.
53604	0.9 g	Single Antibiotic with Neomycin	Bulk	2,000/case	7 lbs.
53310	0.9 g	Single Antibiotic with Bacitracin	144 ct.	12/case	7 lbs.
53300	0.9 g	Single Antibiotic with Bacitracin	Bulk	2,000/case	7 lbs.

Required Minimum Fill for Class A Kits: 10 Uses

Required Minimum Fill for Class B Kits: 25 Uses



Section 6.5 | Antiseptic

Each antiseptic shall meet the requirements of all applicable FDA requirements and shall be contained in an individual-use, non-reusable application containing at least 0.14 fl. oz. (0.5 g) of antiseptic. (Note: Commonly used applicators are swabs and towelettes. Spray containers containing the minimum number of 1/57 oz. (0.5 g) applications required for Class A or Class B kits are acceptable to meet this requirement.)

Order No.	Size	Description	Carton	Case Qty.	Case Wt.
38401	5" x 8"	Antiseptic Wipe	Bulk	2,000/case	17 lbs.
53000	2 oz.	Antiseptic Spray	-	24/case	5 lbs.
34400	5" x 8"	p.a.w.s.® Antimicrobial Wipes	100 ct.	10/case	10 lbs.
34400B	5" x 8"	p.a.w.s.® Antimicrobial Wipes	Bulk	5,000/case	10 lbs.

Required Minimum Fill for Class A Kits: 10 Uses
 Required Minimum Fill for Class B Kits: 50 Uses



Section 6.7 | Burn Treatment

Each burn treatment shall be a water soluble compound packaged in individual-use applications containing at least 1/32 oz. (0.9 g). (Note: Spray containers containing the minimum number of 1/32 oz. (0.9 g) applications required for Class A or Class B kits are acceptable to meet this requirement. Burn treatment, as required here, is intended to address the treatment of minor burns.)

Order No.	Size	Description	Carton	Case Qty.	Case Wt.
50000	0.125 oz.	Burn Gel	25 ct.	24/case	7 lbs.
50000B	0.125 oz.	Burn Gel	Bulk	1,000/case	10 lbs.
50007	0.9 g	Burn Gel	144 ct.	12/case	7 lbs.
50004	2 oz.	Burn Spray	-	24/case	5 lbs.
53404	0.9 g	First Aid & Burn Cream	Bulk	2,000/case	7 lbs.
53405	0.9 g	First Aid & Burn Cream	25 ct.	36/case	4 lbs.
53410	0.9 g	First Aid & Burn Cream	144 ct.	12/case	7 lbs.

Required Minimum Fill for Class A Kits: 10 Uses
 Required Minimum Fill for Class B Kits: 25 Uses



Section 6.14 | Hand Sanitizer

Hand sanitizers shall be water-soluble with a minimum of 62% ethyl alcohol as the active ingredient. Safetec Instant Hand Sanitizer has an active ingredient of 66.5% ethyl alcohol. (Note: Must contain a minimum of at least 1/32 oz. (0.9 g))





Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
17376	0.9 g	Pouch	25 ct.	36/case	4 lbs.
17377	0.9 g	Pouch	144 ct.	12/case	7 lbs.
17378	0.9 g	Pouch	Bulk	2,000/case	7 lbs.
17352	0.125 oz.	Pouch	100 ct.	10/case	12 lbs.
17353	0.057 oz.	Pouch	100 ct.	10/case	7 lbs.

Required Minimum Fill for Class A Kits: 6 Uses
 Required Minimum Fill for Class B Kits: 10 Uses



Is Your Class A First Aid Kit ANSI Compliant?

Class A First Aid Kit Checklist

First Aid Kit Supply	Minimum Quantity	Minimum Size/Vol.	Do You Have It?
Adhesive Bandage	16	1 x 3 in.	
Adhesive Tape	1	2.5 yd (total)	
Antibiotic Application 	10	1/57 oz. (0.5 g)	
Antiseptic 	10	1/57 oz. (0.5 g)	
Breathing Barrier	1	N/A	
Burn Dressing (gel soaked)	1	4 x 4 in.	
Burn Treatment 	10	1/32 oz. (0.9 g)	
Cold Pack	1	4 x 5 in.	
Eye Covering (with means of attachment)	2	2.9 sq. in.	
Eye/Skin Wash	1	1 fl oz. total	
Foil Blanket	1	52 x 84 in.	
First Aid Guide	1	N/A	
Hand Sanitizer 	10	1/32 oz. (0.9 g)	
Medical Exam Gloves	2 pairs	N/A	
Roller Bandage	1	2 in. x 4 yd	
Scissors	1	N/A	
Splint	0	4 x 24 in.	
Sterile Pad	2	3 x 3 in.	
Tourniquet	0	1.5 in. (width)	
Trauma Pad	2	5 x 9 in.	
Triangular Bandage	1	40 x 40 x 56 in.	

*Information based on the ANSI/ISEA Z308.1-2021 American National Standards for Minimum Requirements for Workplace First Aid Kits and Supplies.

 Ready to ship Safetec product



Missing something? We've got you covered.





ANSI compliant, ready to ship Safetec products:

- Triple Antibiotic and Single Antibiotic Ointments
- Antiseptic Wipes and Spray
- First Aid & Burn Cream, Burn Gel, and Burn Spray
- Instant Hand Sanitizer

Note: Safetec does not manufacture pre-packaged first aid kits. However, we can assemble kits using Safetec supplies and/or customer-provided supplies.

Is Your Class B First Aid Kit ANSI Compliant?

Class B First Aid Kit Checklist

First Aid Kit Supply	Minimum Quantity	Minimum Size/Vol.	Do You Have It?
Adhesive Bandage	50	1 x 3 in.	
Adhesive Tape	2	2.5 yd (total)	
Antibiotic Application 	25	1/57 oz. (0.5 g)	
Antiseptic 	50	1/57 oz. (0.5 g)	
Breathing Barrier	1	N/A	
Burn Dressing (gel soaked)	2	4 x 4 in.	
Burn Treatment 	25	1/32 oz. (0.9 g)	
Cold Pack	2	4 x 5 in.	
Eye Covering (with means of attachment)	2	2.9 sq. in.	
Eye/Skin Wash	1	4 fl oz. total	
Foil Blanket	1	52 x 84 in.	
First Aid Guide	1	N/A	
Hand Sanitizer 	20	1/32 oz. (0.9 g)	
Medical Exam Gloves	4 pairs	N/A	
Roller Bandage	2	2 in. x 4 yd	
	1	4 in. x 4 yd	
Scissors	1	N/A	
Splint	1	4 x 24 in.	
Sterile Pad	4	3 x 3 in.	
Tourniquet	1	1.5 in. (width)	
Trauma Pad	4	5 x 9 in.	
Triangular Bandage	2	40 x 40 x 56 in.	

*Information based on the ANSI/ISEA Z308.1-2021 American National Standards for Minimum Requirements for Workplace First Aid Kits and Supplies.

 Ready to ship Safetec product



Need to print more checklists? No worries.

Scan the QR code to download!

Section B2. Each work environment is unique. Accordingly, it is expected that a first aid kit containing the minimum required first aid supplies will be augmented with additional items or additional quantities of required supplies based upon the specific hazards existing in a workplace environment.



Oral Pain Relief

Formulated with Benzocaine for Pain Relief

Safetec's Oral Pain Relief Gel provides rapid and soothing relief from oral discomfort. Formulated with Benzocaine, this peppermint flavored gel effectively numbs and alleviates pain, making it ideal for toothaches, mouth sores, and gum irritation.

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
53117	0.75 g	Pouch	144 ct.	12/case	7 lbs.



1% Hydrocortisone Cream

Temporary Relief of Minor Skin Irritations

Safetec's Hydrocortisone Cream (1%) is designed to temporarily relieve itching associated with minor skin irritations, inflammation, and rashes such as eczema, poison ivy, and dermatitis. Trust it to swiftly address redness, swelling, and itching from various skin irritations, including insect bites.

Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
53104	0.9 g	Pouch	Bulk	2,000/case	7 lbs.
53110	0.9 g	Pouch	144 ct.	12/case	7 lbs.

Sting Relief

with Lidocaine for Fast-Acting Pain Relief

Safetec's Sting Relief Wipes are a must-have for outdoor workers and enthusiasts. Ideal for hiking, camping, or any outdoor activity, these single-use, pre-moistened wipes provide instant pain relief with fast-acting Lidocaine and prevent infection with antiseptic. Individually packaged for portability, they make discomfort from bites and stings disappear with ease.



Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
52001	2 oz.	Spray Bottle	-	24/case	5 lbs.
52013	1.13" x 2.75"	Individual Wipe	Bulk	3,000/case	12 lbs.
52014	1.13" x 2.75"	Individual Wipe	150 ct.	20/case	12 lbs.
52015	1.13" x 2.75"	Individual Wipe	10 ct.	100/case	6 lbs.
52017	1.13" x 2.75"	Individual Wipe	25 ct.	36/case	4 lbs.

52001 is a "Made to Order" only product.

ANSI Compliant First Aid Kit Supplements

SaniZide Pro 1®

1-Minute Surface Disinfectant Spray and Wipes

SaniZide Pro 1® is a broad spectrum, EPA-registered surface disinfectant that combines cleaning and disinfecting in one simple step with its ready-to-use, alcohol-based, hospital-grade formula. SaniZide Pro 1® makes staying compliant even easier by reducing the risk of cross-contamination with true, one-minute kill claims that are effective against 48 microorganisms and safe for repeated use on hard, non-porous surfaces.



Order No.	Size	Product Type	Carton/ Canister	Case Qty.	Case Wt.
35910	32 oz.	Spray Bottle	-	6/case	14 lbs.
35923	6" x 6.75"	Wipe Canister	160 ct.	12/case	22 lbs.
35926	8" x 8"	Individual Wipe	50 ct.	6/case	8 lbs.
35927	8" x 8"	Individual Wipe	Bulk	500/case	15 lbs.
35941	5 Gallon	Carboy	-	1/case	40 lbs.
35955	55 Gallon	Drum	-	1/case	520 lbs.

SaniZide Plus® Germicidal Solution

Kills 99% of Bacteria

SaniZide Plus® is a one-step[†] germicidal cleaner and deodorant designed for general cleaning, disinfecting and controlling non-pathogenic mold and mildew on hard, non-porous inanimate surfaces. Quickly removes dirt, grime, non-pathogenic mold and mildew, food residue, blood and other organic matter commonly found in health care facilities. It also eliminates odors leaving hard non-porous surfaces smelling clean and fresh. Use where odors are a problem. This product is effective against 48 microorganisms on hard non-porous surfaces.

† "One Step" disinfectant claim is limited to association with non-food contact surface use.

* when used according to disinfection directions



Order No.	Size	Product Type	Carton	Case Qty.	Case Wt.
34800	4 oz.	Spray Bottle	-	24/case	8 lbs.
34805	16 oz.	Spray Bottle	-	12/case	15 lbs.
34810	32 oz.	Spray Bottle	-	6/case	15 lbs.
34815	1 Gallon	Refill Bottle	-	4/case	38 lbs.
34826	8" x 8"	Individual Wipe	50 ct.	6/case	13 lbs.
34827	8" x 8"	Individual Wipe	Bulk	1,000/case	32 lbs.
34855	55 Gallon	Drum	-	1/case	520 lbs.

SAMPLE REQUESTS

Safetec's sample program allows you to test any product featured in our catalog. Whether you're stocking your inventory, supporting a customer bid, or exploring private labeling, we're here to help.

How to Order:

- **Contact Your Account Manager:** They'll be happy to submit a request.
- **Visit Our Website** and navigate to the "Distributors" dropdown menu to access the form.
- **Request Via QR Code:** Follow the steps below to submit a sample request now.



Grab your phone & open up the camera



Point it at the QR Code



Tap the notification to open the link

Scan me to request your samples!



While your phone is out, scan below to follow us on social media!







(800) 456-7077



info@safetec.com



www.safetec.com
www.privatelabel.safetec.com



887 Kensington Avenue
Buffalo, NY 14215
United States of America



© 2025 Safetec of America, Inc.
Rev 25.10