

Spill Leader Kit Directions

Kit Number: (15205)

Kit Contents:

- Nitrile Gloves (Quantity: 3 Pairs)
- Red Z[®] Solidifier (5oz. Bottle)
- Green-Z[®] Solidifier (5oz. Bottle)
- Acid Lock[®] Solidifier (5oz. Bottle)
- Alky[®] Solidifier (5oz. Bottle)
- Petro Lock Absorbent (2oz. Bottle)
- Scoop/Scraper
- Zip Lock Bags (Quantity: 6 / Dimensions: 9"x12")
- Hazardous Waste D.O.T. Label (Quantity: 2)
- Non-Hazardous Waste D.O.T. Label (Quantity: 2)

Directions:

1. Determine type of material spilled.
2. Notify proper authorities.
3. Move and keep people away from accident scene.
4. Do not walk into or touch any spilled material.
5. Put on disposable gloves..
6. Place chosen neutralizer and solidifying agent in a circle around the spill.
7. Sprinkle remaining neutralizer and solidifying agent over remaining spill.
8. Remove solidified spill with scoop and scraper(s) and carefully place in the proper bag provided. Disposal bag should be chosen dependent upon spill and your local, state and federal regulations for infected solid waste.
9. Dispose of bag as required by local, state and federal regulations for infected solid waste.