

Instruction Guide

Spill Leader Kit

Reorder # 15205 (Wall Mount)



What's Included:

- Adhesive Wall Mount
- 5 oz. Acid Lock Solidifier
- 5 oz. Alky Solidifier
- 5 oz. Green Z® Solidifier
- 5 oz. Petro Lock Absorbent
- 2 oz. Red Z® Solidifier
- 2 Non-Hazardous Waste D.O.T. Labels
- 6 Scoop/Scrapers
- 6 Zip Lock Bags (9" x 12")
- 2 Hazardous Waste D.O.T. Labels
- 3 Pairs of Nitrile Gloves



Scan here to view the instruction video!

Step 1: Choose Solidifier



Acid Lock

Rapidly neutralizes and solidifies low pH spills.

(Acetic Acid, Hydrochloric Acid, Nitric Acid, and Sulfuric Acid)



Alky

Rapidly neutralizes and solidifies high pH spills.

(Ammonium Hydroxide, Sodium Hydroxide, and Potassium Hydroxide)



Green-Z®

Solidifies blood and body fluids, as well as chemical spills.



Petro Lock


Rapidly absorbs petroleum based spills.

(Gasoline, Solvents, Crude Oil, and Paint Sludge)



Red Z®

Solidifies blood/ body fluids (Contains Chlorine for deodorizing)

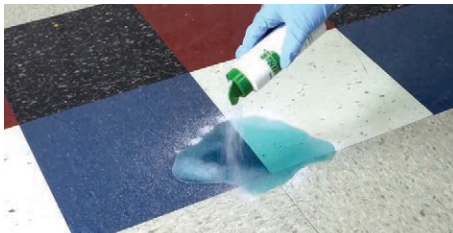

Step 2: Solidifier

Step 2: Sprinkle Solidifier over the spill. Allow liquid to congeal for safer handling and transport.



Step 3: Scoop & Scraper

Step 3: Collect and remove gelled material with Scoop/Scraper.



Step 4: Place gelled material along with Scoop/Scraper into Zip Lock Bag.



Step 5: Wash Hands



Step 6: D.O.T. Label

Step 6: Fill out D.O.T. Label and stick onto Zip Lock Bag. Dispose according to local, state, and federal regulations.

Rev 25.11