

Multi-Purpose Spill Kit



This kit solidifies gallons of various spills. The easy to carry and store molded plastic case fits easily behind a truck seat, or spill response cart. It was developed for use by medical, industrial, transportation, and spill response personnel.

Item#	Description	Case Qty.	Case Wt.
15201	Multi-Purpose Spill Kit	1 each	10 lbs.

Compliance

Meets D.O.T. 1713 hazardous discharge; EPA 263.30 immediate action and 263.31 discharge clean up.

Each Kit Contains:

- Pair of Nitrile Gloves
- Vented Goggles
- Respirator
- Green-Z® Solidifier (15 oz bottle)
- Acid Lock® Solidifier (15 oz bottle)
- Alky® Solidifier (15 oz bottle)
- Petro Lock Absorbent (8 oz bottle)
- Scoop/Scraper (Qty: 3)
- Caution Tape Roll
- Drop Cloth
- Zammie Sewer Stopper (15" x 21")
- Clear Disposal Bags (Qty: 3 – 30" x 37")
- Twist Ties (Qty: 3)
- D.O.T. Incident Report Tag
- Instructions

Spill Leader Kit

Our wall mounted spill kit is designed for emergency response or hazardous lab spills. The solid solutions neutralize, stabilize, and solidify a variety of aqueous spills wherever they occur.



Item#	Description	Case Qty.	Case Wt.
15205	Spill Leader Kit	1 each	5 lbs.

Compliance

Meets D.O.T. 1713 hazardous discharge; EPA 263.30 immediate action and 263.31 discharge clean up.

Each Kit Contains:

- Nitrile Gloves (Qty: 3 pairs)
- Red Z® Solidifier (5 oz bottle)
- Green-Z® Solidifier (5 oz bottle)
- Acid Lock® Solidifier (5 oz bottle)
- Alky® Solidifier (5 oz bottle)
- Petro Lock Absorbent (2 oz bottle)
- Scoop/Scraper (Qty: 6)
- Zip Lock Bags (Qty: 6 – 9" x 12")
- Hazardous Waste D.O.T. Label (Qty: 2)
- Non-Hazardous Waste D.O.T. Label (Qty: 2)

See other side for solidifier descriptions

Spill Cleanup Solidifiers

Safely handle a variety of spills within your facility using these Safetec Solidifiers



Red Z® Solidifier (contains chlorine)

Solidifies potentially infectious blood, and body fluids

Sprinkle Red Z® evenly over the spill. Allow sufficient time for solidification. Remove solidified material with scoop and scraper. Clean and disinfect contaminated area after use and wash thoroughly.*



Green-Z® Solidifier

Solidifies potentially infectious blood, body fluids and chemical spills

Sprinkle Green-Z® to surround and evenly distribute over the spill. Allow sufficient time for solidification. Remove solidified material with scoop and scraper. Clean and disinfect contaminated area after use and wash thoroughly.*

**Always wear protective clothing when cleaning up and/or handling a spill. Dispose of infectious and hazardous waste in accordance with all local, state and federal regulations.*

To watch a demonstration on how to use these products visit safetec.com/resources/



Safetec of America, Inc.

887 Kensington Avenue Buffalo, NY 14215
1-800-456-7077 • www.safetec.com



Alky® Solidifier

Rapidly neutralizes and solidifies high pH spills

- Ammonium hydroxide
- Sodium Hydroxide
- Potassium hydroxide

Sprinkle Alky® directly onto the spill, covering it completely. The spill will turn blue, then yellow when it has safely neutralized. Once it becomes a semi-solid gel, it is ready for cleanup with a scoop and scraper.*



Acid Lock® Solidifier

Rapidly neutralizes and solidifies low pH spills

- Acetic acid
- Hydrochloric acid
- Nitric acid
- Sulfuric acid

Sprinkle Acid Lock® directly onto the spill, covering it completely. The spill will turn pink, then yellow when it has safely neutralized. Once it becomes a semi-solid gel, it is ready for cleanup with a scoop and scraper.*



Petro Lock Solidifier

Rapidly absorbs petroleum based spills

- Gasoline
- Solvents
- Crude oil
- Paint sludge

Sprinkle Petro Lock directly onto the spill, covering it completely to easily absorb fluid. Once it becomes a semi-solid gel, it is ready for cleanup with a scoop and scraper.*

(Petro Lock is oil specific and will not absorb water base liquids)